

Lovibond® Water Testing

Tintometer® Group



Step 1

Please open the instruction „HexLoad_de_gb.pdf“ in the „HexLoad“ dir.
Print the instruction and have it at hand when updating.

Step 2

To update your SpectroDirect, MultiDirect, PoolDirect, PC9in1, TurbiDirect, PCPECTRO II or MaxiDirect open the program HEXLoad.

Step 3

Additional every Product has its own firmware. You will find it in the folder „Firmware“ fitting to your product.

Step 4

Update your device as described in the instruction „HexLoad_de_gb.pdf“.

Step 5

Download the current version of the instruction for your product from our website.

Instruction Software-Update

Lovibond® Water Testing

Tintometer® Group



Changelog

New regarding MultiDirect, MaxiDirect, PoolDirect & PC9in1

- new Cyanuric Acid Method (CyA-TEST) replaces Cyanuric Acid Tablet
Important Note: After the Update, the Cyanuric Acid Tablet can not be used any more. For more Information, visit our Website.
 - OTZ Function (OneTimeZero)
 - New Methods MaxiDirect
 - Alkalinity-M HR with tablet reagents
 - Chloride with liquid reagent
 - Chlorine HR with powder reagent
 - Chlorine HR DPD with tablet reagents
 - Iron with liquid reagent
 - Molybdate LR with powder reagent / liquid reagent
 - Phosphate HR with liquid reagent
 - Polyacrylate with liquid reagent
 - Triazole with powder reagent
 - New Methods MultiDirect
 - Alkalinity-M HR with tablet reagents
 - Chlorine HR DPD with tablet reagents
 - Molybdate LR with powder reagent / liquid reagent
-

New regarding SpectroDirect

- new Cyanuric Acid Method (CyA-TEST) replaces Cyanuric Acid Tablet
Important Note: After the Update, the Cyanuric Acid Tablet can not be used any more. For more Information, visit our Website.
- New Methods SpectroDirect
 - Molybdate LR with powder reagent / liquid reagent
 - Monochloramin
 - Phosphonate
 - Suspended Solids
 - Alkalinity-M HR with tablet reagents
 - Chlorine HR DPD with tablet reagents